

Denver Instrument
DI352 Titrator
(800) 321-1135 x 262
Jim Rinehart

Vitamin C 0.031804% ± 0.00041

Printed the 6/5/2006
3:53:08PM


Fruit Salad Juice

| | | | |
|----------------|---------------------|----------------|--------------|
| Operator | Jim Rinehart | Method ID | VIT C |
| Workstation | DI352 | Method type | Titration |
| Coupled method | No | Titration type | Monotonic IP |
| Analysis date | 6/5/2006 11:57:13AM | Electrode ID | DI300746.1 |
| Status | OK | Titrant | I2 0.1 |
| | | Tests | 3 |

Sample

| | | | |
|-----------|------------------|----------------|------|
| Sample ID | Dole Fruit Salad | Advised amount | 30 g |
| Type | Normal | | |
| Batch no. | None | | |

Curves

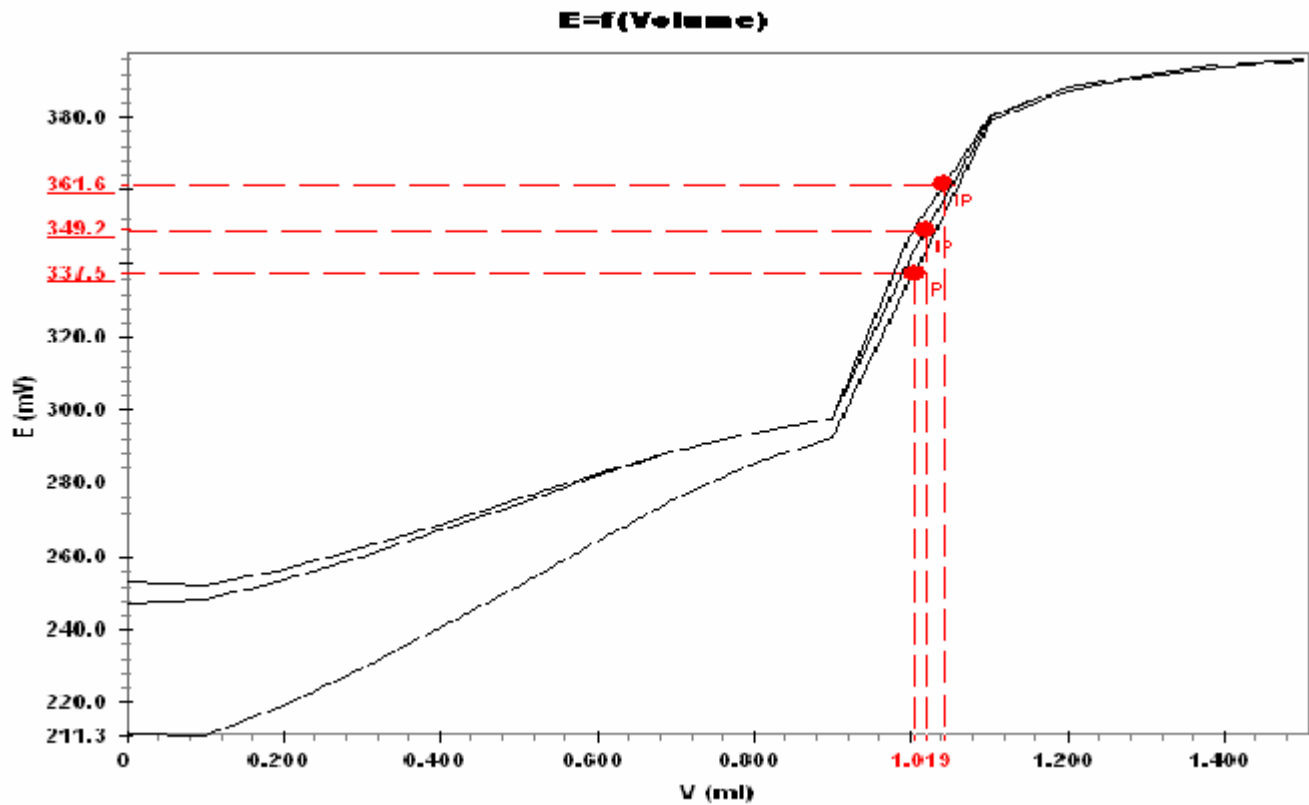

Result

| Name | Value | Tests | Standard dev | Acceptation | T°C |
|---------------|------------|-------|--------------|-------------|------|
| Ascorbic Acid | 0.031804 % | 3 | 0.0004118 | OK | 25.0 |

Comments**Signature**

